

E-Line 80 Refractometer



User Guide



**Bellingham
+ Stanley**

E-Line 80 Refractometer User Guide (Eng)

B + S Code : 44-893

Issue 1C

January 2010

Table I

<i>Order Code</i>	<i>Range</i>	<i>Scale Division</i>	<i>Scale No.</i>	<i>Scale Ranges</i>
44-807	E-Line 80	0.5 °Brix minimum	1	0-50 °Brix
			2	50-80 °Brix

© Copyright Bellingham + Stanley Ltd 2010

Every effort has been made to ensure the accuracy of the contents of this manual. However, Bellingham + Stanley Ltd can assume no responsibility for errors contained in the manual or their consequences.

Bellingham + Stanley pursues a policy of continuous product development and improvement and as such, information given in this manual may be updated or withdrawn without notice. Some publicity materials prior to January 2010 will not reflect such product improvements, noticeably the change in colour of the instrument barrel from black to bright silver. This change does not affect the performance of the instrument under normal operating conditions.

Printed in United Kingdom.

Bellingham+Stanley Ltd.

Longfield Road, Tunbridge Wells,
Kent, TN2 3EY, United Kingdom
Phone: +44 (0) 1892 500400
Fax: +44 (0) 1892 543115
sales@bellinghamandstanley.co.uk

Bellingham+Stanley Inc.


1000 Hurricane Shoals Road, Building D,
Suite 300, Lawrenceville, GA 30043, USA
Phone: 770 822 6898
Fax: 770 822 9165
sales@bs-rfm-inc.com

Product Description

The E-Line 80 is a dual scale Brix refractometer ideal for use in the food industry and other applications requiring concentration measurements over a wide measurement range.

Applying the sample to the refractometer:

Lift the illuminating prism¹ and apply the sample to the measurement prism². Close the prism and view the borderline (light/dark demarcation line) through the eyepiece³.

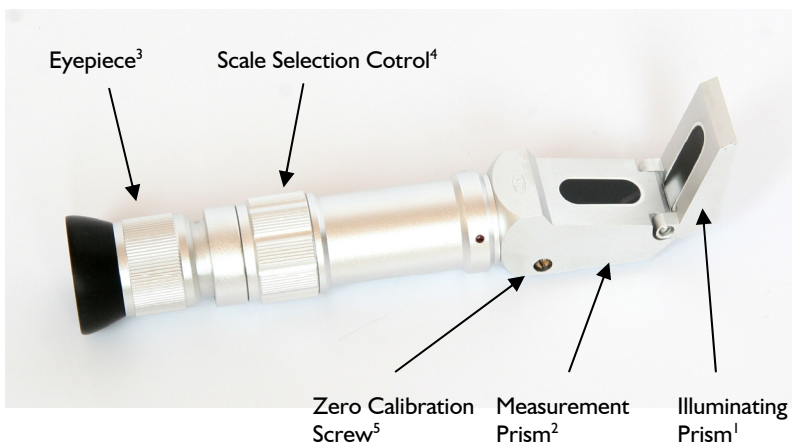

Optical glass is relatively soft and care should be taken not to scratch the prism surface. Do not use metal spatulas or glass rods to apply samples but instead use softer materials such as plastic.

Always check sample Health & Safety Data before applying to the refractometer.

When applying samples to the prism which are likely to cause harm to skin or eyes, wear appropriate protective clothing and glasses.

Taking a reading:

Depending on the concentration of the sample, select the range of measurement^(Table 1) by adjusting the scale selection control⁴. If the concentration is unknown, apply the sample and step through the ranges until a clear borderline is found.

The concentration of the sample can be measured at the point the borderline crosses the scale. Take care to ensure that the reading is taken from the correct scale range.

Refractive Index values can be obtained by using a conversion table available from Bellingham + Stanley Ltd, see Table 3.

Cleaning the prism:

Thoroughly clean the prism after use with water or other suitable solvent and dry with clean tissue. The prism surface could be damaged by strong alkalis or acids if left in contact for long periods of time. Clean samples from the prism as soon as practicable. Wiping the prism surface occasionally with alcohol will remove any build-up of oils left from the samples.

Correcting the reading for temperature:

The refractometer is calibrated to read correctly at 20°C. The refractive index of the measured sample will vary with temperature and the reading must be corrected to take account of this. Please use the Table 2 below to adjust your reading accordingly.

Temperature Correction Table (Table 2)

This table gives mass fraction corrections to refractometric tables for sucrose solutions at 589nm for temperature different from 20°C. Taken from ICUMSA, Appendix 2, SPS-3 (1998) page 8.

Temp. (°C)	Measured Sucrose Brix (Mass Fraction)																	
	0	5	10	15	20	25	30	35	40	45	50	55	60	65	70	75	80	85
15	-0.29	-0.30	-0.32	-0.33	-0.34	-0.35	-0.36	-0.37	-0.37	-0.38	-0.38	-0.38	-0.38	-0.38	-0.38	-0.38	-0.37	-0.37
16	-0.24	-0.25	-0.26	-0.27	-0.28	-0.28	-0.29	-0.30	-0.30	-0.30	-0.31	-0.31	-0.31	-0.31	-0.31	-0.30	-0.30	-0.30
17	-0.18	-0.19	-0.20	-0.20	-0.21	-0.21	-0.22	-0.22	-0.23	-0.23	-0.23	-0.23	-0.23	-0.23	-0.23	-0.23	-0.23	-0.22
18	-0.12	-0.13	-0.13	-0.14	-0.14	-0.14	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15	-0.15
19	-0.06	-0.06	-0.07	-0.07	-0.07	-0.07	-0.07	-0.08	-0.08	-0.08	-0.08	-0.08	-0.08	-0.08	-0.08	-0.08	-0.08	-0.07
20	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00	0.00
21	0.06	0.07	0.07	0.07	0.07	0.07	0.08	0.08	0.08	0.08	0.08	0.08	0.08	0.08	0.08	0.08	0.08	0.07
22	0.13	0.14	0.14	0.14	0.15	0.15	0.15	0.15	0.16	0.16	0.16	0.16	0.16	0.16	0.15	0.15	0.15	0.15
23	0.20	0.21	0.21	0.22	0.22	0.23	0.23	0.23	0.23	0.24	0.24	0.24	0.24	0.23	0.23	0.23	0.23	0.22
24	0.27	0.28	0.29	0.29	0.30	0.30	0.31	0.31	0.31	0.32	0.32	0.32	0.32	0.31	0.31	0.31	0.30	0.30
25	0.34	0.35	0.36	0.37	0.38	0.38	0.39	0.39	0.40	0.40	0.40	0.40	0.40	0.39	0.39	0.38	0.38	0.37
26	0.42	0.43	0.44	0.45	0.46	0.46	0.47	0.47	0.48	0.48	0.48	0.48	0.48	0.47	0.47	0.46	0.46	0.45
27	0.50	0.51	0.52	0.53	0.54	0.55	0.55	0.56	0.56	0.56	0.56	0.56	0.56	0.55	0.55	0.54	0.53	0.52
28	0.58	0.59	0.60	0.61	0.62	0.63	0.64	0.64	0.64	0.65	0.65	0.64	0.64	0.63	0.63	0.62	0.61	0.60
29	0.66	0.67	0.68	0.70	0.71	0.71	0.72	0.73	0.73	0.73	0.73	0.73	0.72	0.72	0.71	0.70	0.69	0.67
30	0.74	0.76	0.77	0.78	0.79	0.80	0.81	0.81	0.82	0.82	0.81	0.81	0.80	0.80	0.79	0.78	0.76	0.75
31	0.83	0.84	0.85	0.87	0.88	0.89	0.89	0.90	0.90	0.90	0.90	0.89	0.89	0.88	0.87	0.86	0.84	0.82
32	0.92	0.93	0.94	0.96	0.97	0.98	0.98	0.99	0.99	0.99	0.99	0.98	0.97	0.96	0.95	0.93	0.92	0.90
33	1.01	1.02	1.03	1.05	1.06	1.07	1.07	1.08	1.08	1.08	1.07	1.07	1.06	1.04	1.03	1.01	1.00	0.98
34	1.10	1.11	1.13	1.14	1.15	1.16	1.16	1.17	1.17	1.16	1.16	1.15	1.14	1.13	1.11	1.09	1.07	1.05
35	1.19	1.21	1.22	1.23	1.24	1.25	1.25	1.26	1.26	1.25	1.25	1.24	1.23	1.21	1.19	1.17	1.15	1.13
36	1.29	1.30	1.31	1.33	1.34	1.34	1.35	1.35	1.35	1.34	1.34	1.33	1.31	1.29	1.28	1.25	1.23	1.20
37	1.39	1.40	1.41	1.42	1.43	1.44	1.44	1.44	1.44	1.43	1.43	1.41	1.40	1.38	1.36	1.33	1.31	1.28
38	1.49	1.50	1.51	1.52	1.53	1.53	1.54	1.54	1.53	1.53	1.52	1.50	1.48	1.46	1.44	1.42	1.39	1.36
39	1.59	1.60	1.61	1.62	1.63	1.63	1.63	1.63	1.63	1.62	1.61	1.59	1.57	1.55	1.52	1.50	1.47	1.43
40	1.69	1.70	1.71	1.72	1.73	1.73	1.73	1.73	1.72	1.71	1.70	1.68	1.66	1.63	1.61	1.58	1.54	1.51

Calibration:

The E-line 80 is calibrated to read accurately at 20°C. Readings may be taken at any other temperature by applying a correction factor (Table 2).

Alternatively the E-line 80 has the facility to correct for temperature by manual offset. Prior to taking a reading, ensure the temperature of the instrument is stable. Apply distilled water and check that the reading is zero (Scale 1). If the reading is not zero, use the screwdriver supplied to adjust the zero calibration screw⁵ so that the borderline crosses the zero mark on the scale. No temperature correction is required for this method but as the relationship between refractive index and temperature is not linear, small errors may occur at the top end of the scale compared to readings near water.

Care Warning:



This refractometer is a precision optical instrument and should be handled with care. Do not drop or subject it to sharp knocks.

°Brix to Refractive Index Conversion Table (Table 3)

The table below can be used to convert °Brix readings taken from the instrument scale into refractive index.

°Brix	Refractive Index at 589.3nm and 20.0°C	°Brix	Refractive Index at 589.3nm and 20.0°C	°Brix	Refractive Index at 589.3nm and 20.0°C
0	1.33299	30	1.38115	60	1.44193
1	1.33442	31	1.38296	61	1.44420
2	1.33586	32	1.38478	62	1.44650
3	1.33732	33	1.38661	63	1.44881
4	1.33879	34	1.38846	64	1.45113
5	1.34026	35	1.39032	65	1.45348
6	1.34175	36	1.39220	66	1.45584
7	1.34325	37	1.39409	67	1.45822
8	1.34477	38	1.39600	68	1.46061
9	1.34629	39	1.39792	69	1.46303
10	1.34782	40	1.39986	70	1.46546
11	1.34937	41	1.40181	71	1.46790
12	1.35093	42	1.40378	72	1.47037
13	1.35250	43	1.40576	73	1.47285
14	1.35408	44	1.40776	74	1.47535
15	1.35568	45	1.40978	75	1.47787
16	1.35729	46	1.41181	76	1.48040
17	1.35891	47	1.41385	77	1.48295
18	1.36054	48	1.41592	78	1.48552
19	1.36218	49	1.41799	79	1.48811
20	1.36384	50	1.42009	80	1.49071
21	1.36551	51	1.42220		
22	1.36720	52	1.42432		
23	1.36889	53	1.42647		
24	1.37060	54	1.42862		
25	1.37233	55	1.43080		
26	1.37406	56	1.43299		
27	1.37582	57	1.43520		
28	1.37758	58	1.43743		
29	1.37936	59	1.43967		

Data Source: ICUMSA Methods Book, Specification and Standard SPS-3 (2000), Refractometry and Tables - (Official)

Page intentionally left blank

Page intentionally left blank